

# **Human Chorionic Gonadotropin (hCG) Enzyme Immunoassay Test**

Cat. No: KBBA09

Ver2.0

Immunoassay for the quantification of hCG in human Serum

*For Research Purposes Only. Not for use in diagnostic or therapeutic procedures. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited.*

**Introduction:**

Human chorionic gonadotropin (hCG) is a glycoprotein hormone produced by the placenta. During normal pregnancy the level of this hormone in serum and urine gradually increases up to about the eighth week of pregnancy. hCG has two subunits, namely  $\alpha$  and  $\beta$ . The  $\alpha$ -subunit is similar to the  $\alpha$ -subunit of the anterior pituitary gland glycoprotein hormones, namely, LH, TSH and FSH. In normal pregnancy the increase in hCG starts at about the 5th day after conception and continues to rise until it reaches a maximum at about the eighth week. It is a well known fact that hCG is also a tumour marker which is very important in the diagnosis of choriocarcinoma.

**Intended Use:**

This kit is intended for determination of Human Chorionic Gonadotropin Hormone (hCG) in human serum.

**Principle:**

The method employs the quantitative sandwich enzyme immunoassay technique. Antibodies to hCG are pre-coated onto microwells. Samples and standards are pipetted into microwells and anti hCG antibodies linked to HRP is added to the wells. hCG present in the sample and standards will bind to the plate and HRP Conjugate simultaneously. Washing is done to remove the unbound samples, standards and HRP conjugate. The ready to use substrate solution (TMB) is added to microwells and color develops directly proportional to the amount of hCG in the standards and sample. Color development is then stopped by addition of stop solution. Absorbance is measured at 450 nm.

Please note that a two-step procedure is included and is to be used for assaying serum from pregnant women.

**Materials Provided:**

1. Anti-hCG coated Microtiter plate (96 wells) – 1 no
2. Control 1 – 0.5 ml
3. Control 2 - 0.5ml
4. hCG Standard 0 IU/L – 2ml
5. hCG Standard (0.5 ml/vial) - 2.5, 10, 25, 100 and 500 IU/L
6. Anti-hCG Antibody:HRP Conjugate – 300  $\mu$ l/vial
7. Assay Buffer – 25 ml
8. Wash Buffer – 50 ml
9. TMB Substrate – 16 ml
10. Stop Solution – 6 ml
11. Instruction Manual

**Materials to be provided by the End-User:**

1. Microtiter Plate Reader able to measure absorbance at 450 nm.
2. Adjustable pipettes and multichannel pipettor to measure volumes ranging from 25 $\mu$ l to 1000 $\mu$ l
3. Deionised (DI) water
4. ELISA Plate shaker
5. Wash bottle or automated microplate washer
6. Log-Log graph paper or software for data analysis

**Handling/Storage:**

1. All reagents should be stored at 2<sup>o</sup>C to 8<sup>o</sup>C for stability.
2. All the reagents and wash solutions are stable until the expiration date of the kit.
3. Before using, bring all components to Room Temperature (18-25 °C). Upon assay completion return all components to appropriate storage conditions.
4. The Substrate is light-sensitive and should be protected from direct sunlight or UV sources.

**Health Hazard Warnings:**

1. This kit contains human material. The source material used for manufacture of this kit tested negative for HBsAg, HIV 1/2 and HCV by FDA-approved methods. However, no method can completely assure absence of these agents. Therefore, all human blood products, including serum samples, should be considered potentially infectious. Handling should be as defined by an appropriate national biohazard safety guideline or regulation, where it exists
2. Reagents that contain preservatives may be harmful if ingested, inhaled or absorbed through the skin.
3. For Research or Manufacturing use only.

**Reagent Preparation (all reagents should be diluted immediately prior to use):**

1. Label any aliquots made with the kit Lot No and Expiration date and store it at appropriate conditions mentioned.
2. Bring all reagents to Room Temperature before use.
3. To make Wash Buffer (1X); dilute 50ml of 10X Wash Buffer in 450 ml of DI water.
4. Anti-hCG Antibody:HRP Conjugate; dilute 240µl of Anti-hCG antibody: HRP in 12ml of Assay Buffer.

**Specimen Collection and Preparation**

Serum should be prepared from a whole blood specimen obtained by acceptable medical techniques. This kit is for use with serum samples without additives only.

**Assay Procedure:****One-Step Procedure:**

1. Bring all reagents to room temperature prior to use. It is strongly recommended that all Standards and Samples be run in duplicates or triplicates. A standard curve is required for each assay. All steps must be performed at room temperature (RT).
2. Pipette out 25µl of **Standards, Controls** and **Samples** into the respective wells as mentioned in the work list.
3. Pipette 100µl of the anti-hCG-HRP Conjugate into each well.
4. Incubate on a plate shaker (approximately 200 rpm) for 1 hour at room temperature.
5. Aspirate and wash plate 3 times with **Wash Buffer (1X)** and blot residual buffer by firmly tapping plate upside down on absorbent paper.
6. Add 100µl of **TMB Substrate** in to each well.
7. Incubate the plate on a plate shaker for 10-15 minutes in dark at RT. **DO NOT SHAKE** or else it may result in higher backgrounds and worse precision. Positive wells should turn bluish in color.
8. Pipette out 50µl of **Stop Solution**. Wells should turn from blue to yellow in color.
9. Read the absorbance at 450 nm with a microplate reader within 20 minutes after addition of the stop solution.

**Two Step Procedure (Note: this procedure is to be used when assaying serum from pregnant women)**

1. Bring all reagents to room temperature prior to use. It is strongly recommended that all Standards and Samples be run in duplicates or triplicates. A standard curve is required for each assay. All steps must be performed at room temperature (RT).
2. Pipette out 25µl of **Standards, Controls** and **Samples** into the respective wells as mentioned in the work list.
3. Pipette 100µl of assay buffer into each well.
4. Incubate on a plate shaker (approximately 200 rpm) for 1 hour at room temperature.
5. Aspirate and wash plate 3 times with **Wash Buffer (1X)** and blot residual buffer by firmly tapping plate upside down on absorbent paper.
6. Pipette 100µl of the anti-hCG-HRP conjugate into each well.
7. Incubate on a plate shaker (approximately 200 rpm) for 1 hour at room temperature.
8. Aspirate and wash plate 3 times with **Wash Buffer (1X)** and blot residual buffer by firmly tapping plate upside down on absorbent paper.
9. Pipette 100µl of TMB substrate into each well.
10. Incubate on a plate shaker for 10-15 minutes at room temperature.
11. Pipette 50µl of stopping solution into each well.
12. Read the plate on a microwell plate reader at 415nm within 20 minutes after addition of the stopping solution.

**Example of a Work list**

<b>Well #</b>	<b>Contents</b>	<b>Abs at 450nm</b>	<b>Mean Absorbance</b>	<b>IU/L hCG equiv.</b>
1A	Zero Std			
2A	Zero Std			
1B	2.5IU/L			
2B	2.5IU/L			
1C	10IU/L			
2C	10IU/L			
1D	25IU/L			
2D	25IU/L			
1E	100IU/L			
2E	100IU/L			
1F	500IU/L			
2F	500IU/L			
1G	Control 1			
2G	Control 1			
1H	Control 2			
2H	Control 2			

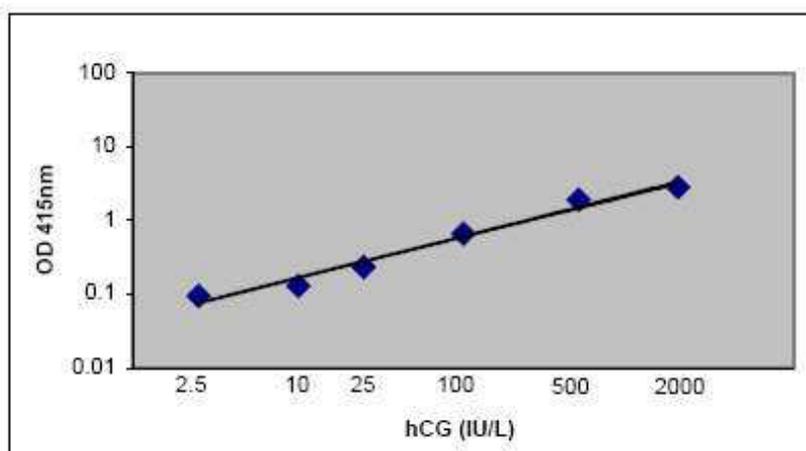
**Calculation of Results:**

Determine the Mean Absorbance for each set of duplicate or triplicate Standards, Controls and Samples. Subtract the Mean Absorbance of the Zero Standards (background) from each well. Using log-log graph paper or a computer program, plot the optical densities of each standard on the Y-axis versus the corresponding concentration of the standards on the X-axis. Draw the best fit straight line through the standard points. Draw the best fit straight line through the standard points. To determine the unknown hCG concentrations, find the unknown's Mean Absorbance value on the Y-axis and draw a horizontal line to the standard curve. At the point of intersection, draw a vertical line to the X-axis and read the hCG Concentration. Computer based curve-fitting software may be preferred.

If a sample reads more than 500 IU/L then dilute it with Standard 0 at a dilution of no more than 1:20. The result should be multiplied with the dilution factor.

**Typical Data:**

This standard curve was generated at KRISHGEN for demonstration purposes only. A standard curve must be run with each assay.

**Performance Characteristics:****1. Sensitivity:**

- Limit Of Detection:** It is defined as the lowest detectable concentration corresponding to a signal of Mean of '0' standard plus  $2 \times SD$ . 10 replicates of '0' standards were evaluated and the LOD was found to be 0.7 IU/L.
- Limit of Quantitation:** It is defined as the lowest concentration for which Coefficient of Variation is  $<20\%$ . The LOQ is found to be  $<2.5$  IU/L.

**Precautions:**

Do not mix reagents from different kits or lots. Reagents and/or antibodies from different manufacturers should not be used with this set.

**Quality Control:**

It is recommended that for each laboratory assay appropriate quality control samples in each run to be used to ensure that all reagents and procedures are correct.

**Troubleshooting**

<b>Problem</b>	<b>Possible cause</b>	<b>Investigation/Actions</b>
High Absorbances	1. Cross-contamination from other specimens	> Repeat assay taking care when washing and pipetting.
	2. Insufficient or inefficient washing or reading	> Check washer efficiency
	3. Wavelength of filter not correct.	> Check that the wavelength is 450nm. If a dual wavelength spectrophotometer is available, set the reference filter between 600-650 nm.
	4. High assay background.	> Repeat assay and include a well that contains only sample diluent or sample absorbent (i.e. a blank well).
	5. Contaminated TMB	> Check that TMB is colorless or faint blue.
	6. Incubation time too long or incubation temperature too high.	> Check incubation time and temperature. > Check incubator is at the correct temperature.
	7. Incorrect dilution of serum	> Repeat assay, ensuring correct serum dilution is used.
Low Absorbances	1. Incubation time too short or incubation temperature too low.	> Ensure time and temperature of assay incubation are correct
	2. Incorrect dilution or pipetting of sera	> Check incubator is set at the correct temperature.
	3. Incorrect filter wavelength.	> Repeat assay ensuring correct dilutions and volumes are used
	4. Contaminated Conjugate solution.	> Ensure controls are sufficiently mixed. > Check the wavelength is set at 450nm. If a dual wavelength spectrophotometer is available, set the reference filter between 600-650nm. > Dispense conjugate directly from the bottle using clean pipette tip; avoid transferring Conjugate to another container if possible. > Do not return unused Conjugate to bottle. > Ensure all pipettes and probes used to dispense the Conjugates are clean and free from serum, detergent and bleach.
	5. Kit has expired.	> Check expiration date of kit and do not use if expired.
	6. Air blank reading high	> Investigate causes of high background absorbance.
	7. Incorrect storage of kit.	> Ensure kit is stored at 2-8°C, plate is sealed in foil pouch and desiccant sachet is blue/purple.
	8. Kit reagents not equilibrated at room temperature	> Allow sufficient time for reagents to equilibrate to room temperature prior to assay.
	9. Incorrect reagents used.	> Check the reagents used match those listed on the specification sheet.
	10. Over washing of plate (e.g. inclusion of a long soak step).	> Repeat assay using recommended wash procedure.
Poor Duplicates	1. Poor mixing of samples.	> Mix reagents gently and equilibrate to room temperature.
	2. Poor pipette precision	> Calibration may need to be checked. > Check pipetting technique-change pipette tip for each sample and ensure excess liquid is wiped from the outside of the tip.
	3. Addition of reagents at inconstant timing intervals; reagent addition takes too long, air bubbles when adding reagents.	> Use consistent timing when adding reagents. > Ensure all dilutions are made before commencing addition to plate.
	4. Inefficient washing - Wash buffer left in wells, inconsistent washing, inadequate washing.	> Improve pipetting technique and skill. > Tap out wash buffer after washing. > Check wells are sufficiently and uniformly filled and aspirated when washing.
	5. Reader not calibrated or warmed up prior to plate reading.	> Check reader precision > Check reader manual to ascertain warm up time of instrument.
	6. Optical pathway not clean	> Gently wipe bottom of plate. > Check reader light source and detector are clean.
	7. Spillage of liquid from wells	> Repeat assay, taking care not to knock the plate or splash liquid
	8. Serum samples exhibit microbial growth, haemolysis or lipaemia.	> It is not recommended to use serum samples exhibiting microbial
	9. Uneven well volumes due to evaporation.	> Cover plate with a lid or plate sealer (not provided).

- |                    |   |   |
|--------------------|---|---|
| All wells Yellow   | <ol style="list-style-type: none"> <li>1. Contaminated TMB.</li> <li>2. Contaminated reagents (e.g. Conjugate, Wash buffer)</li> <li>3. Incorrect dilution of serum.</li> <li>4. Incorrect storage of kit.</li> <li>5. Inefficient washing- Wash buffer left in wells, inconsistent washing, inadequate washing.</li> <li>6. If conjugate reconstitution is required- Conjugate reconstituted incorrectly.</li> </ol> | <ul style="list-style-type: none"> <li>&gt; Check TMB is colorless or faint blue.</li> <li>&gt; Check reagents for turbidity.</li> <li>&gt; Repeat assay, ensuring correct serum dilution is used.</li> <li>&gt; Ensure kit is stored at 2-8°C, plate is sealed in foil pouch and desiccant sachet is blue / purple.</li> <li>&gt; Tap out wash buffer after washing.</li> <li>&gt; Check wells are sufficiently and uniformly filled an aspirated washing.</li> <li>&gt; Repeat assay ensuring Conjugate is reconstituted according to assay method.</li> </ul>  |
| All wells Negative | <ol style="list-style-type: none"> <li>1. Test not performed correctly – correct reagents not added or not added in the correct sequence</li> <li>2. Contaminated Conjugate solution.</li> <li>3. Over- washing of plate (e.g. inclusion of a long soak step).</li> <li>4. Incorrect storage of kit.</li> <li>5. Wash Buffer made up with Stop Solution instead of Wash Buffer Concentrate</li> </ol>                 | <ul style="list-style-type: none"> <li>&gt; Check procedure and check for unused reagents.</li> <li>&gt; Ensure that Stop Solution was not added before Conjugate or TMB.</li> <li>&gt; Ensure that serum was diluted in correct Sample diluent; e.g. do not use Sample Absorbent for an IgG ELISA.</li> <li>&gt; Dispense Conjugate directly from the bottle using a clean pipette tip; avoid transferring Conjugate to another container if possible.</li> <li>&gt; Do not return unused Conjugate to bottle.</li> <li>&gt; Ensure all pipettes and probes used to dispense the Conjugate are clean and free from serum, detergent and bleach.</li> <li>&gt; Repeat assay using recommended wash procedure.</li> <li>&gt; Ensure kit is stored at 2-8°C, plate is sealed in foil pouch and desiccant sachet is blue / purple.</li> <li>&gt; Ensure Wash Buffer is made up correctly.</li> </ul> |

**LIMITED WARRANTY**

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any) from the maker thereof of such non Krishgen made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen Biosystems shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

*Krishgen Biosystems. 2014*

**THANK YOU FOR USING KRISHGEN PRODUCT!**