

# KRIBIOLISA™ Denosumab (PROLIA™) ELISA

**REF** : KBI1026

Ver 2.1

**RUO**

Enzyme Immunoassay for the quantitative determination Denosumab (PROLIA™) in serum, plasma and cell culture supernatant

<b>RUO</b>	For Research Use	<b>REF</b>	Catalog Number
	Store At	<b>LOT</b>	Batch Code
	Manufactured By		Biological Risk
	Expiry Date		Consult Operating Instructions

*For Research Use Only. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited.*

**REF** KBI1026

 96 tests



**KRISHGEN BioSystems** | For US / Europe: toll free +1(888)-970-0827 tel: +1(562)-568-5005

For Asia / India: tel: +91(22)-49198700

Email: sales@krishgen.com

**Introduction:**


Denosumab (trade names Prolia and Xgeva) is a human monoclonal antibody for the treatment of osteoporosis, treatment-induced bone loss, metastases to bone, and giant cell tumor of bone. Denosumab is contraindicated in people with low blood calcium levels. The most common side effects are joint and muscle pain in the arms or legs. Denosumab is a RANKL inhibitor, which works by preventing the development of osteoclasts which are cells that break down bone (bone resorption).

**Intended Use:**

The KRIBIOLISA™ Denosumab (PROLIA™) ELISA is used as an analytical tool for quantitative determination of Denosumab (PROLIA™) in serum, plasma and cell culture supernatant.

**Principle:**

The method employs the quantitative sandwich enzyme immunoassay technique. Antibodies to Denosumab are pre-coated onto microwells. Samples and standards are pipetted into microwells and human Denosumab present in the sample are bound by the capture antibody. Then, a HRP (horseradish peroxidase) conjugated Anti-Denosumab antibody is pipetted and incubated. After washing microwells in order to remove any non-specific binding, the ready to use substrate solution (TMB) is added to microwells and color develops proportionally to the amount of Denosumab in the sample. Color development is then stopped by addition of stop solution. Absorbance is measured at 450 nm.

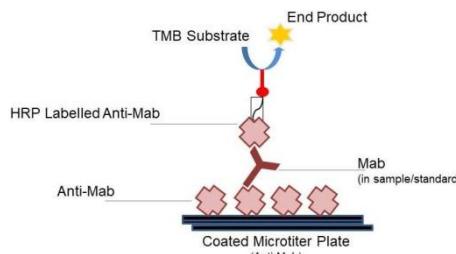**Materials Provided:**

Part	Description	Qty
Anti-Denosumab Coated Microtiter Plate	96 well polystyrene microplate (12 strips of 8 wells) coated with Anti-Denosumab monoclonal antibody.	1 x 96 wells
Denosumab Standard	Recombinant Denosumab in a buffered protein base with preservative sodium azide <0.1% (lyophilized, 1 ug/ml)	2 vials
Anti-Denosumab:HRP Conjugate	Anti-Denosumab conjugated to Horseradish Peroxidase with protein stabilizer and preservatives 0.02% methylisothiazolone and 0.02% bromonitrodioxane.	12 ml
(1X) Sample Diluent	Buffered protein base with preservative sodium azide <0.1%	50 ml
(1X) Standard Diluent	Buffered protein base with 1:10 dilution human serum and preservative sodium azide < 0.1%	10 ml
(20X) Wash Buffer	20-fold concentrated solution of buffered surfactant with preservative thiomersol < 0.01%. May turn yellow over time.	25 ml
TMB Substrate	Stabilized Chromogen	12 ml
Stop Solution	0.73M Phosphoric Acid	12 ml
Instruction Manual		1 no

**Materials to be provided by the End-User:**

1. Microtiter Plate Reader able to measure absorbance at 450 nm.
2. Adjustable pipettes and multichannel pipettor to measure volumes ranging from 25 ul to 1000 ul
3. Deionized (DI) water
4. Wash bottle or automated microplate washer
5. Graph paper or software for data analysis
6. Timer
7. Absorbent Paper

**Handling/Storage:**

1. All reagents should be stored at 2°C to 8°C for stability.
2. All the reagents and wash solutions should be used within 12 months from manufacturing date.
3. Before using, bring all components to room temperature (18-25°C). Upon assay completion ensure all components of the kit are returned to appropriate storage conditions.
4. The Substrate is light-sensitive and should be protected from direct sunlight or UV sources.

**Health Hazard Warnings:**

1. Reagents that contain preservatives may be harmful if ingested, inhaled or absorbed through the skin.
2. For Research Use Only.

**Sample Preparation and Storage:**

Blood is taken by venipuncture. Serum is separated after clotting by centrifugation. Plasma can be used, too. Lipaemic, hemolytic or contaminated samples should not be run. Repeated freezing and thawing should be avoided. If samples are to be used for several assays, initially aliquot samples and keep at - 20°C.

For Cell Culture Supernatant – If necessary, centrifuge to remove debris prior to analysis. Samples can be stored at -20°C or -80°C. Avoid repeated freeze-thaw cycles.

**Preparation Before Use:**

Allow samples to reach room temperature prior to assay. Take care to agitate patient samples gently in order to ensure homogeneity.

Test Sample preparation - Samples have to be diluted 1:10 (v/v), e.g. 1 ul sample + 9 ul Sample Diluent prior to assay. The samples may be kept at 2 - 8°C for up to three days. Long-term storage requires -20°C.

**Reagent Preparation (all reagents should be diluted immediately prior to use):**

1. Label any aliquots made with the kit Lot No and Expiration date and store it at appropriate conditions mentioned.
2. Bring all reagents to Room temperature before use.
3. To make Wash Buffer (1X); dilute 25 ml of 20X Wash Buffer in 475 ml of DI water.
4. **Standards Preparation:** Reconstitute the lyophilized standard in 1 ml Standard diluent to get a concentration of 1 ug/ml. Dilute 160 ul of original **Standard (1 ug/ml)** with 840 ul of Standard Diluent to generate a **160 ng/ml Standard Solution**. Prepare further **Standards** by diluting the Standard Solution as per the below table. Use the Standard Diluent as the Zero Standard (Standard No.0).

Standard Concentration	Standard Vial	Dilution Particulars
1 ug/ml	Lyophilized Standard	Lyophilized standard + Reconstitute in 1 ml Standard Diluent
160 ng/ml	Standard No.5	160 ul Reconstituted Standard (1 ug/ml) + 840 ul Standard Diluent
80 ng/ml	Standard No.4	500 ul Standard No.5 + 500 ul Standard Diluent
40 ng/ml	Standard No.3	500 ul Standard No.4 + 500 ul Standard Diluent
20 ng/ml	Standard No.2	500 ul Standard No.3 + 500 ul Standard Diluent
10 ng/ml	Standard No.1	500 ul Standard No.2 + 500 ul Standard Diluent
0 ng/ml	Standard No.0	Only Standard Diluent

Use the Standards immediately upon reconstitution. Discard balance standard after use. Do not store them for further experiments

**Procedural Notes:**

1. In order to achieve good assay reproducibility and sensitivity, proper washing of the plates to remove excess un-reacted reagents is essential.

2. High Dose Hook Effect may be observed in samples with very high concentrations of Denosumab (PROLIA). High Dose Hook Effect is due to excess of antibody for very high concentrations of Denosumab (PROLIA) present in the sample. High Dose Hook effect is most likely encountered from samples early in the purification process. If Hook Effect is possible, the samples to be assayed should be diluted with a compatible diluent. Thus if the Denosumab (PROLIA) concentration of the undiluted sample is less than the diluted sample, this may be indicative of the Hook Effect.
3. Avoid assay of Samples containing sodium azide ( $\text{NaN}_3$ ), as it could destroy the HRP activity resulting in under-estimation of the amount of Denosumab (PROLIA).
4. It is recommended that all Standards and Samples be assayed in duplicates.
5. Maintain a repetitive timing sequence from well to well for all the steps to ensure that the incubation timings are same for each well.
6. If the Substrate has a distinct blue color prior to use it may have been contaminated and use of such substrate can lead to compromisation of the sensitivity of the assay.
7. The plates should be read within 30 minutes after adding the Stop Solution.
8. Make a work list in order to identify the location of Standards and Samples.

#### Assay Procedure:

1. It is strongly recommended that all Controls and Samples be run in duplicates or triplicates. A standard curve is required for each assay. All steps must be performed at 37°C
2. Pipette **100 ul** of **Standards** or **Samples** into the respective wells.
3. Cover the plate and incubate for 60 minutes at 37°C
4. Aspirate and wash plate 4 times with **Wash Buffer (1X)** and blot residual buffer by firmly tapping plate upside down on absorbent paper. Wipe of any liquid from the bottom outside of the microtiter wells as any residue can interfere in the reading step.
5. Pipette without delay in the same order **100 ul** of **Anti-Denosumab:HRP Conjugate** into each well.
6. Cover the plate and incubate for 60 minutes at 37°C
7. Aspirate and wash plate 4 times with **Wash Buffer (1X)** and blot residual buffer by firmly tapping plate upside down on absorbent paper. Wipe of any liquid from the bottom outside of the microtiter wells as any residue can interfere in the reading step.
8. Add **100 ul** of **TMB Substrate** in each well.
9. Incubate the plate at 37°C for 30 minutes in dark. DO NOT SHAKE or else it may result in higher backgrounds and worse precision. Positive wells should turn bluish in color.
10. Pipette out **100 ul** of **Stop Solution**. Wells should turn from blue to yellow in color.
11. Read the absorbance at 450 nm with a microplate reader.

#### Calculation of Results:

Determine the Mean Absorbance for each set of duplicate or triplicate Standards and Samples. Using Semi-Log graph paper, plot the average value (absorbance 450nm) of each standard on the Y-axis versus the corresponding concentration of the standards on the X-axis. Draw the best fit curve through the standard points. To determine the unknown Denosumab concentrations, find the unknown's Mean Absorbance value on the Y-axis and draw a horizontal line to the standard curve. At the point of intersection, draw a vertical line to the X-axis and read the Denosumab Concentration. If samples were diluted, multiply by the appropriate dilution factor. Software which is able to generate a cubic spline curve-fit is best recommended for automated results.


#### Note:

It is recommended to repeat the assay at a different dilution factor in the following cases:

- If the sample absorbance value is below the first standard.
- If the absorbance value is equivalent or higher than the 160 ng/ml standard.

Standard Concentration (ng/ml)	Abs A	Abs B	Mean Abs	Interpolated Concentration	% Interpolated Concentration against Actual Concentration
0	0.057	0.059	0.058	--	--
10	0.276	0.284	0.280	10.6	106.4
20	0.500	0.572	0.536	20.7	103.7
40	0.911	1.056	0.984	38.4	95.9
80	1.846	1.912	1.879	81.4	101.7
160	3.121	2.566	2.843	159.4	99.6

Typical Graph

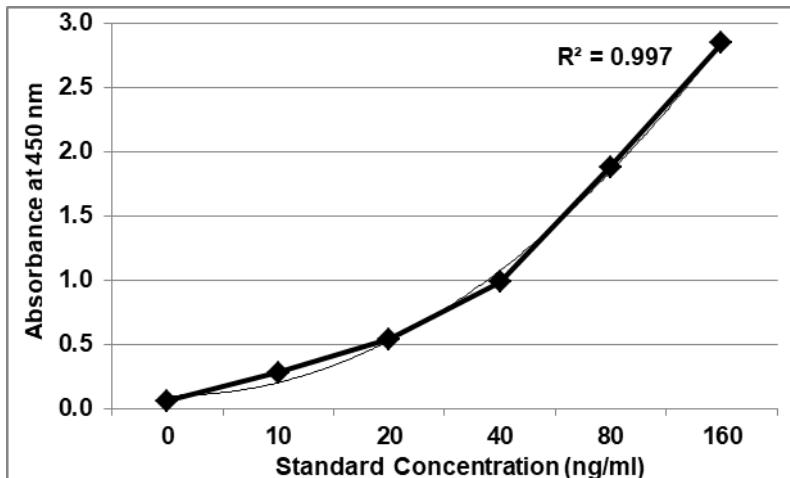

#### Quality Control:

It is recommended that for each laboratory assay appropriate quality control samples in each run to be used to ensure that all reagents and procedures are correct.

#### Performance Characteristics of the Kit:

This kit has been validated as per EMA/FDA guidelines in line with ICH Code for Harmonization of Biological Assays.

#### Sensitivity:

**Limit Of Detection:** It is defined as the lowest detectable concentration corresponding to a signal of Mean of '0' standard plus 2\* SD.

10 replicates of '0' standards were evaluated and the LOD was found to be 5 ng/ml

#### Linearity:

Standards provided in the kit will be used for measuring the linearity range of Denosumab present in matrix. The standard has been calibrated and validated against commercially sourced Prolia Injection<sup>1</sup>.

#### Precision:

Precision is defined as the percent coefficient of variation (%CV) i.e. standard deviation divided by the mean and multiplied by 100. Assay precision was determined by both intra (n=5 assays) and inter assay (n=5 assays) reproducibility on two pools with low (10ng/ml), medium (40ng/ml) and high (160ng/ml) concentrations. While actual precision may vary from laboratory to laboratory and technician to technician, it is recommended that all operators achieve precision below these design goals before reporting results.

Pool	Intra Assay %CV	Inter Assay %CV
Low	<12%	<15%
Medium	<10%	<12%
High	<10%	<12%

- **This kit is For Research Use only.** Follow the working instructions carefully.
- The expiration dates stated on the kit are to be observed. The same relates to the stability stated for reagents
- Do not use or mix reagents from different lots.
- Do not use reagents from other manufacturers.
- Avoid time shift during pipetting of reagents.
- All reagents should be kept in the original shipping container.
- Some of the reagents contain small amount of sodium azide (< 0.1 % w/w) as preservative. They must not be swallowed or allowed to come into contact with skin or mucosa.
- Source materials maybe derived from human body fluids or organs used in the preparation of this kit were tested and found negative for HBsAg and HIV as well as for HCV antibodies. However, no known test guarantees the absence of such viral agents. Therefore, handle all components and all patient samples as if potentially hazardous.
- Since the kit contains potentially hazardous materials, the following precautions should be observed
  - Do not smoke, eat or drink while handling kit material
  - Always use protective gloves
  - Never pipette material by mouth
  - Wipe up spills promptly, washing the affected surface thoroughly with a decontaminant.
- In any case GLP should be applied with all general and individual regulations to the use of this kit.

<sup>1</sup>Prolia Injection is the registered trade mark of Amgen Inc.



## SCHEMATIC ASSAY PROCEDURE

1. Remove all components, 30 minutes before adding into the assay plate.



2. Avoid repeated cool-thaw of the components as there will be a loss of activity and this can affect the results.



3. Pipette **100 µl Standards / Samples** into each well.

4. Cover plate and incubate for at 37°C.

5. Aspirate and wash wells 4 times with **Wash Buffer (1X)**.

6. Pipette **100 µl Anti-Denosumab:HRP** into each well.

7. Cover plate and incubate for at 37°C.

8. Aspirate and wash wells 4 times with **Wash Buffer (1X)**.

9. Pipette **100 µl TMB Substrate** into each well.

10. Cover plate and incubate for at 37°C.

11. Pipette **100 µl Stop Solution** into each well.

12. Read absorbance at 450nm with a microplate reader within of stopping reaction.

## Typical Example of a Work List

Well #	Contents	Absorbance at 450nm	Mean Absorbance	ng/ml Denosumab (PROLIA) equivalent
1A	zero std			
2A	zero std			
1B	10 ng/ml			
2B	10 ng/ml			
1C	20 ng/ml			
2C	20 ng/ml			
1D	40 ng/ml			
2D	40 ng/ml			
1E	80 ng/ml			
2E	80 ng/ml			
1F	160 ng/ml			
2F	160 ng/ml			
1G	Sample			
2G				
1H	Sample			
2H				
3A	Sample			
4A				
3B	Sample			
4B				

**LIMITED WARRANTY**

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any) from the maker thereof of such non Krishgen made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

**THIS WARRANTY IS EXCLUSIVE.** The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen Biosystems shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, and strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

Krishgen Biosystems. 2023

**THANK YOU FOR USING KRISHGEN PRODUCT!**

KRISHGEN BIOSYSTEMS®, GENLISA®, DHARMAPLEX™, GENBULK™, GENLISA™, KRISHZYME®, KRISHGEN®, KRIBIOLISA®, KRISHPLEX®, TITANIUM®, QUALICHEK® are registered trademarks of KRISHGEN BIOSYSTEMS. ©KRISHGEN BIOSYSTEMS. ALL RIGHTS RESERVED.

**KRISHGEN BIOSYSTEMS | OUR REAGENTS | YOUR RESEARCH |**

## SYMBOLS KEY

	<b>Anti-Denosumab Coated Microtiter Plate (12x8 wells)</b>
	<b>Denosumab Standard , lyophilized</b>
	<b>Conjugate Horseradish Peroxidase</b>
	<b>(1X) Sample Diluent</b>
	<b>(1X) Standard Diluent</b>
	<b>(20X) Wash Buffer</b>
	<b>TMB Substrate</b>
	<b>Stop Solution</b>
	<b>Consult Instructions for Use</b>
	<b>Catalog Number</b>
	<b>Expiration Date</b>
	<b>Storage Temperature</b>