






GENLISA™ Human Anti-SARS-CoV-2 (Covid-19) IgG nucleocapsid protein Qualitative ELISA

REF : KBVH015-3

Ver 2.1


RUO

Enzyme Immunoassay for the Qualitative Estimation of IgG Antibodies to Human SARS-CoV-2 (Covid-19) in human serum and plasma, respiratory specimens and cell culture supernatant

RUO	For Research Use Only	REF	Catalog Number
	Store At	LOT	Batch Code
	Manufactured By		Biological Risk
	Expiry Date		Consult Operating Instructions

For Research Use Only. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited.

REF KBVH015-3

 96 tests

KRISHGEN BioSystems For US/Europe Customers: toll free +1(888)-970-0827 | tel +1(562)-568-5005
For Asia/India Customers: tel +91(22)-49198700
Email: sales@krishgen.com | http://www.krishgen.com

Introduction:

The GENLISA™ ELISA kits are used for assessing the specific biomarker in samples analytes which may be serum, plasma and cell culture supernatant as validated with the kit. The kit employs a sandwich ELISA technique which leads to a higher specificity and increased sensitivity compared to conventional competitive ELISA kits which employ only one antibody.

There is no standard reference SARS-CoV-2 antigen material available; accordingly, absolute analytical sensitivity cannot be calculated. The kit uses a polyclonal antibody as Positive Control. The assay allows the qualitative determination of samples.

Intended Use:

The GENLISA™ Human Anti-SARS-CoV-2 (Covid-19) IgG ELISA kit is used as an analytical tool for qualitative estimation of IgG antibodies to Human SARS-CoV-2 (Covid-19) in human serum and plasma, respiratory specimens and cell culture supernatant.

Principle:

The method employs sandwich ELISA technique. Human SARS-CoV-2 protein is pre-coated onto microwells. Samples and controls are pipetted into microwells and IgG Antibodies to human SARS-CoV-2 (Covid-19) present in the sample are bound by the protein antigen. After incubation the wells are washed and followed by HRP-conjugated Detection IgG Antibody is pipetted and incubated to form a complex. After washing microwells in order to remove any non-specific binding, the substrate solution (TMB) is added to the microwells. Color develops proportionally to the amount of IgG Anti-Human SARS-CoV-2 (Covid-19) in the samples. Color development is then stopped by addition of stop solution. Absorbance is measured at 450 nm.

Materials Provided:

1. Recombinant SARS-CoV-2 (Covid-19) nucleocapsid protein Coated Microtiter Plate (12 x 8 wells) - 1 no
2. Positive Control - 0.5 ml vial
3. Negative Control - 0.5 ml vial
4. Anti-Human IgG:HRP Conjugate - 12 ml
5. (5X) Assay Diluent - 25 ml
6. (20X) Wash Buffer - 25 ml
7. TMB Substrate - 12 ml
8. Stop Solution - 12 ml

Materials to be provided by the End-User:

1. Microtiter Plate Reader able to measure absorbance at 450 nm.
2. Adjustable pipettes and multichannel pipettor to measure volumes ranging from 25 ul to 1000 ul
3. Deionized (DI) water
4. Wash bottle or automated microplate washer
5. Graph paper or software for data analysis
6. Timer
7. Absorbent Paper

Handling/Storage:

1. Store main kit components at 2-8°C.
2. Before using, bring all components to room temperature (18-25°C). Upon assay completion return all components to appropriate storage conditions.
3. The Substrate is light-sensitive and should be protected from direct sunlight or UV sources.

Health Hazard Warnings:

Reagents that contain preservatives may be harmful if ingested, inhaled or absorbed through the skin.

**Sample Preparation and Storage:**

Specimens should be clear and non-hemolyzed. Samples should be run at a number of dilutions to ensure accurate quantitation.

Blood is taken by venipuncture. Serum is separated after clotting by centrifugation. Plasma can be used, too. Lipaemic, hemolytic or contaminated samples should not be run. Repeated freezing and thawing should be avoided.

Samples should be diluted 1:5000 (v/v) for optimal recovery, (for example 1 ul sample + 4999 ul sample diluent) prior to assay. In cases where matrix interferences is under or over observed, the samples may be diluted with Sample Diluent accordingly.

The samples may be kept at 2 - 8°C for up to three days. For long-term storage please store at -20°C.

Note: *Grossly hemolyzed samples are not suitable for use in this assay.*

Cell Culture Supernates - Remove particulates by centrifugation and assay immediately or aliquot and store samples at -20°C or lower temperature. Avoid repeated freeze-thaw cycles. If the use of original supernate sample or low dilution (<5 fold) are necessary due to the expected low concentration of antigen supernates need be adjust to neutral pH condition before assay.

Should you desire to inactivate the virus, use a (5X) Lysis Buffer (optional, not provided). Add 1/5 volume of (5X) Lysis Buffer to sample (i.e. add 50 ul (5X) Lysis Buffer to 200 ul sample). Vortex well.

Note:

The sample should be diluted to within the working range of the assay in 1X Assay Diluent. The exact dilution must be determined based on the concentration of specific target in individual samples.

Respiratory Sample - Centrifuge samples for 20 minutes at 10000 x g at 2-8°C. Collect supernatant and carry out the assay immediately.

Reagent Preparation (all reagents should be diluted immediately prior to use):

1. Label any aliquots made with the kit Lot No and Expiration date and store it at appropriate conditions mentioned.
2. Bring all reagents to Room Temperature before use.
3. To make **Assay Diluent (1X)**; dilute **25 ml of 5X Assay Diluent in 100 ml of DI water**.
4. To make **Wash Buffer (1X)**; dilute **25 ml of 20X Wash Buffer in 475 ml of DI water**.

Procedural Notes:

1. In order to achieve good assay reproducibility and sensitivity, proper washing of the plates to remove excess un-reacted reagents is essential.
2. Avoid assay of Samples containing sodium azide (NaN_3), as it could destroy the HRP activity resulting in erroneous results for the presence of Anti-Human SARS-CoV-2 (Covid-19).
3. It is recommended that the Controls and Samples be assayed in duplicates.
4. Maintain a repetitive timing sequence from well to well for all the steps to ensure that the incubation timings are same for each well.
5. If the Substrate has a distinct blue color prior to use it may have been contaminated and use of such substrate can lead to compromisation of the sensitivity of the assay.
6. The plates should be read within 30 minutes after adding the Stop Solution.
7. Make a work list in order to identify the location of Controls and Samples.

Assay Procedure:

1. Pipette 100 ul of **Controls** and **diluted Samples** to the respective wells. Seal plate and incubate for 1 hour at Room Temperature (18-25°C).
2. Aspirate and wash plate 4 times with **Wash Buffer (1X)** and blot residual buffer by firmly tapping plate upside down on absorbent paper. Wipe of any liquid from the bottom outside of the microtiter wells as any residue can interfere in the reading step. All the washes should be performed similarly.
3. Add 100 ul of **Anti-Human IgG:HRP Conjugate** to each well.
4. Seal plate and incubate for 1 hour at Room Temperature (18-25°C).
5. Wash plate 4 times with **Wash Buffer (1X)** as in step 2.
6. Pipette 100 ul of **TMB Substrate solution**.
7. Incubate in the dark for 15 minutes at Room Temperature.
8. Stop reaction by adding 100 ul of **Stop Solution** to each well.
9. Read absorbance at 450 nm within 30 minutes of stopping reaction.

Interpretation of Results:

Calculation for Cut off Values

Read the sample and negative control wells on microtitre plate reader at 450nm. The OD (Optical Density) of NC (Negative Control) in duplicate should be used for calculating the mean and standard deviation. This is the $Negative_{mean}$.

The cut-off for Positives is equal to a value greater than ($Negative_{mean} + 3*Standard\ Deviation$).

Formula:

Positive Sample Value = OD > ($Negative_{mean} + SD$)

Typical example –

Sample Type	Absorbance #1	Absorbance #2	Mean
Negative	0.131	0.128	0.129
Standard Deviation	$0.131-0.129$ = 0.002	$0.128-0.129$ = -0.001	

$$Mean\ Standard\ Deviation = \sqrt{(0.002)^2 + (-0.001)^2} / 2 = 0.0014$$

$$\begin{aligned} \text{Therefore Cut-off} &= \text{Mean} + 3*SD \\ &= 0.129 + 3* 0.0014 \\ &= 0.129 + 0.0042 \\ &= 0.133 \end{aligned}$$

Positive Samples	> Cut Off
Negative Samples	<= Cut Off
Positive Control Value	> Cut Off
Negative Control Value	Absorbance < 0.5

Quality Control:

It is recommended that for each laboratory assay appropriate quality control samples in each run to be used to ensure that all reagents and procedures are correct. The test is only for screening purpose only and the results should not be the sole basis for clinical diagnosis and treatment. The confirmation of infection with novel coronavirus (COVID-19) must be combined with the patient's clinical symptoms in conjunction to other tests.

In the first week of the onset of the infection with the novel coronavirus (COVID-19) patients results may be negative for IgG. In addition, patients with low immunity or other diseases that affect immune function, failure of important systemic organs, and use of drugs that suppress immune function can also lead to negative results of new coronavirus IgG. Patients previously infected with SARS or other coronavirus strain may cause a light IgG positive in view of similarity of different strains.

Bacterial or fungal contamination of serum specimens or reagents, or cross-contamination between reagents may cause erroneous results.

Performance Characteristics of the Kit:

Sensitivity:

Limit Of Detection: There is no standard reference SARS-CoV-2 antigen material available; accordingly, absolute analytical sensitivity cannot be calculated. The sensitivity adopted using the KBVH015-9 Quantitative ELISA kit was 2 AU/ml. (for more details, please refer the IFU/Kit Insert for KBVH015-9).

Specificity:

Reactivity/Inclusivity

Mutations in the SARS-CoV-2 genome have been identified as the virus has spread, but no serologically unique strains have been described relative to the originally isolated virus (this research is limited at present).

Cross-reactivity of non-SARS-CoV-2 specific antibodies against SARS-CoV-2 nucleocapsid proteins in Anti-SARS-CoV-2 ELISA KBVH015-3 was examined using sera with known antibodies against confirmed past infections.

N	Antibody Positive Sera	Anti-SARS-CoV-2 ELISA #KBVH015-3
1	Beta Corona HKU1*	Negative
4	VCV	Negative
5	HCV	Negative
4	HAV	Negative
3	HBV	Negative
4	EBV	Negative
5	CMV	Negative
5	HSV	Negative

*The patient was tested PCR positive for Beta Corona HKU1 and PCR negative for SARS-CoV-2. Four weeks after PCR testing a serum sample was drawn from the patient and found to be negative in the Anti-SARS-CoV-2 ELISA.

Traceability:

There are no reference standards for SARS-Cov-2 Antibody. The results are reported as positive or negative based on the cut-offs calculated.

Positive Control:

Positive Control provided in the kit was calibrated using the KBVH015-9 ELISA kit. The Value is 15 AU/ml.

Precision:

Precision is defined as the percent coefficient of variation (%CV) i.e. standard deviation divided by the mean and multiplied by 100. Assay precision was determined by both intra (n=5 assays) and inter assay (n=5 assays) reproducibility on two pools with low (3.75 AU/ml), medium (15 AU/ml) and high (30 AU/ml) concentrations.

While actual precision may vary from laboratory to laboratory and technician to technician, it is recommended that all operators achieve precision below these design goals before reporting results.

Pool	Intra Assay %CV	Inter Assay %CV
Low	<10%	<10%
Medium	<5%	<5%
High	<5%	<5%

Recovery:

Human sera and plasma were measured with two replicates and two runs (n = 5). The human sera and plasma were pooled patient and single donor spiked samples. Samples were measured using one lot of reagent. All data met our acceptance criteria for % CV and 95% (CI) Confidence Intervals for % CV.

Matrix	Recovery Range %
Serum (n=5)	87 - 112
Plasma EDTA (n=5)	85 - 114
Plasma Heparin (n=5)	86 - 114

Note: Serum and Plasma were diluted using Standard Diluent provided with the kit

Assay Comparison:

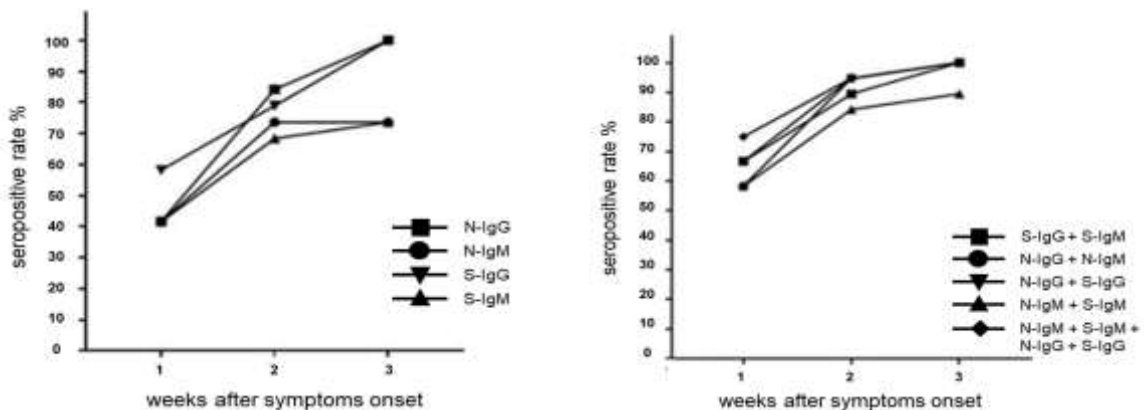
To demonstrate the identity of the Anti-SARS-CoV-2 ELISA and a commercially available COVID-19 IgG / IgM Rapid Test, 9 serum samples were measured using both assays. In comparison, the results of the IgG tests show an increased sensitivity of the anti-SARS-CoV-2 ELISA.

Sample No	Anti-SARS-CoV-2 ELISA	COVID-19 IgG / IgM Rapid Test
1	positive +	negative
2	positive +	positive
3	positive +	negative
4	negative -	negative
5	borderline +	slightly positive
6	positive +++	positive
7	negative -	negative
8	negative -	negative
9	positive ++	positive

Longitudinal Reponse of Antibodies to SARS-CoV-2.

Antibodies response to the Nucleocapsid Proteins and the Spike Proteins of IgG and IgM:

figure a and b



Weeks	N-IgM	N-IgG	S-IgM	S-IgG	N-IgM + N-IgG	S-IgM + S-IgG	N-IgM + S-IgM	N-IgG + S-IgG	N-IgM+ S-IgM+ N-IgG+ S-IgG
1	41.7	41.7	41.7	58.3	58.3	66.7	58.3	66.7	75
2	73.7	84.2	68.4	78.9	94.7	89.5	84.2	94.7	94.7
3	73.7	100	73.7	100	100	100	89.5	100	100

Reference:

EMERGING MICROBES AND INFECTIONS JOURNAL

Accepted author version posted online: 01 May 2020

Kinetics of SARS-CoV-2 specific IgM and IgG responses in COVID-19 patients

Baoqing Sun, Ying Feng, Xiaoneng Mo, Peiyan Zheng, Qian Wang, Pingchao Li, Ping Peng, Xiaoqing Liu, Zhilong Chen, Huimin Huang, Fan Zhang, Wenting Luo, Xuefeng Niu, Peiyu Hu, Longyu Wang, Hui Peng, Zhifeng Huang, Liqiang Feng, Feng Li, Fuchun Zhang, Fang Li, Nanshan Zhong & Ling Chen

<https://doi.org/10.1080/22221751.2020.1762515>

Safety Precautions:

- **This kit is For Research Use Only.** Follow the working instructions carefully.
- The expiration dates stated on the kit are to be observed. The same relates to the stability stated for reagents
- Do not use or mix reagents from different lots.
- Do not use reagents from other manufacturers.
- Avoid time shift during pipetting of reagents.
- All reagents should be kept in the original shipping container.
- Some of the reagents contain small amount of sodium azide (< 0.1 % w/w) as preservative. They must not be swallowed or allowed to come into contact with skin or mucosa.
- Source materials maybe derived from human body fluids or organs used in the preparation of this kit were tested and found negative for HBsAg and HIV as well as for HCV antibodies.
- However, no known test guarantees the absence of such viral agents. Therefore, handle all components and all patient samples as if potentially hazardous.
- Since the kit contains potentially hazardous materials, the following precautions should be observed
 - Do not smoke, eat or drink while handling kit material
 - Always use protective gloves
 - Never pipette material by mouth
 - Wipe up spills promptly, washing the affected surface thoroughly with a decontaminant.
- In any case GLP should be applied with all general and individual regulations to the use of this kit.



References:

Structure, function, and antigenicity of the SARS-CoV-2 nucleocapsid glycoprotein

AC Walls, YJ Park, MA Tortorici, A Wall, AT McGuire...

- Cell, 2020 - Elsevier

Preliminary identification of potential vaccine targets for the COVID-19 coronavirus (SARS-CoV-2) based on SARS-CoV immunological studies

SF Ahmed, AA Quadeer, MR McKay

- Viruses, 2020 - mdpi.com

Interaction between heptad repeat 1 and 2 regions in nucleocapsid protein of SARS-associated coronavirus: implications for virus fusogenic mechanism and identification of ...

S Liu, G Xiao, Y Chen, Y He, J Niu, CR Escalante...

- The Lancet, 2004 - Elsevier

Temporal profiles of viral load in posterior oropharyngeal saliva samples and serum antibody responses during infection by SARS-CoV-2: an observational cohort ...

KKW To, OTY Tsang, WS Leung, AR Tam...

- The Lancet Infectious ..., 2020 - Elsevier

Modulation of the unfolded protein response by the severe acute respiratory syndrome coronavirus nucleocapsid protein

CP Chan, KL Siu, KT Chin, KY Yuen, B Zheng...

- Journal of ..., 2006 - Am Soc Microbiol

Functional assessment of cell entry and receptor usage for SARS-CoV-2 and other lineage B betacoronaviruses

M Letko, A Marzi, V Munster

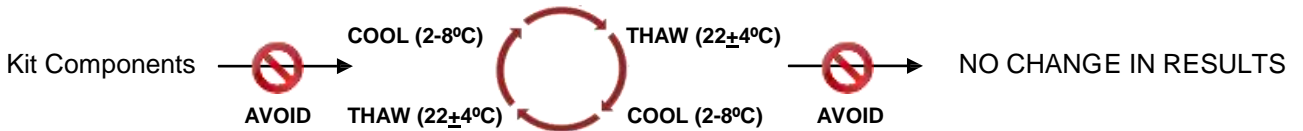
- Nature microbiology, 2020 - nature.com

SCHEMATIC ASSAY PROCEDURE

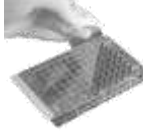

1. Remove all components, 30 minutes before adding into the assay plate.





2. Avoid repeated cool-thaw of the components as there will be a loss of activity and this can affect the results.




3.  Pipette **100 ul Controls and diluted Samples** into the respective wells.

4.  Cover plate and **incubate** for  at Room Temperature.

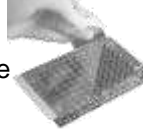

5.  Aspirate and wash wells 4 times with **Wash Buffer (1X)**.

6.  Pipette **100 ul Anti-Human IgG:HRP Conjugate** into each well.

7.  Cover plate and **incubate** for  at Room Temperature.

8.  Aspirate and wash wells 4 times with **(1X) Wash Buffer**.

9.  Pipette **100 ul TMB Substrate** into each well.

10.  Cover plate and **incubate** for  at Room Temperature.

11.  Pipette **100 ul Stop Solution** into each well.

12. Read absorbance at 450nm with a  microplate reader within  of stopping reaction.

Typical Example of a Work List

Well #	Contents	Absorbance at 450nm	Mean Absorbance	Results
1A 2A	+ve Control +ve Control			
1B 2B	-ve Control -ve Control			
1C 2C	<i>Sample</i>			
1D 2D	<i>Sample</i>			
1E 2E	<i>Sample</i>			
1F 2F	<i>Sample</i>			
1G 2G	<i>Sample</i>			
1H 2H	<i>Sample</i>			
3A 4A	<i>Sample</i>			
3B 4B	<i>Sample</i>			

LIMITED WARRANTY

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any) from the maker thereof of such non Krishgen made products or components.

This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen Biosystems shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, and strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

Krishgen Biosystems. 2020

THANK YOU FOR USING KRISHGEN PRODUCT!