

# GENLISA™ Mouse Mycoplasma Pulmonis ELISA

Cat. No: KLM300  
Ver1.3

Enzyme Immunoassay for qualitative screening of antibodies  
against Mouse Mycoplasma Pulmonis in mouse sera

	<b>For Research Use Only</b>		<b>Catalog Number</b>
	<b>Store At</b>		<b>Batch Code</b>
	<b>Manufactured By</b>		<b>Biological Risk</b>
	<b>Expiry Date</b>		<b>Consult Operating Instructions</b>

*For Use On Research Animals Only. Not for use in diagnostic or therapeutic procedures. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited*

**KRISHGEN BioSystems** | Unit Nos3318/319, Shah & Nahar, Off Dr E Moses Road, Worli, Mumbai 400 018.  
Tel: (022) 49198 700 Email: sales@krishgen.com

**Introduction:**

Mycoplasma is a bacteria lacking cell wall. Mycoplasma pulmonis is a species specific organism carried by nearly all pet rats and wild house mice. It colonizes the luminal surface of the respiratory system. It is responsible for respiratory and genital infections in pet rats. It is non-pathogenic to humans and even though it can be carried by humans in nasal passages for short period of time, you yourself cannot be infected. In addition, they are very fragile organisms and rarely remain viable outside the host for long periods making it nearly impossible for it to be transmitted through fomites (e.g., equipments in cages, or carried on clothing). However, how rapidly transmission occurs may vary among rats within a colony.

**Intended Use:**

The Mouse Mycoplasma pulmonis GENLISA™ ELISA is used as an analytical tool for qualitative laboratory screening of presence or absence of antibodies against mouse Mycoplasma pulmonis in the serum of mouse.

**Principle:**

The micro test wells are coated in alternating columns with Mycoplasma pulmonis (MPUL) Antigen and Control-Antigen. During the first incubation with the diluted specimen, any antibodies that are reactive with the *MPUL* Antigen or Control-Antigen will bind to the coated wells. Wash to remove the rest of the sample. Further addition of Enzyme Conjugate it will bind to antibodies which has bound to the antigen coated wells. Unbound HRP conjugate will get removed by washing step after incubation. Then addition of Substrate develops blue color during incubation period and reaction will get stop after addition of stop solution with development of yellow color. The color development is proportional to the concentration of antibodies against mouse *MPUL* in the sample.

**Materials Provided:**

1. Microtitre coated plate (96 wells) - 1 no  
2 holders containing each of 12 positive and 12 negative coated strips
2. Positive Control - 1 ml
3. Negative Control - 1 ml
4. Enzyme Conjugate - 11 ml
5. (20X) Wash Buffer - 2 x 25 ml
6. Sample Diluent - 1 x 30 ml
7. TMB Substrate - 6 ml
8. Stop Solution - 6 ml
9. Instruction Manual

**Materials to be provided by the End-User:**

1. Microplate Reader able to measure absorbance at 405 nm.
2. Adjustable pipettes to measure volumes ranging from 50µl to 1000µl.
3. Deionized (DI) water.
4. Wash bottle or automated microplate washer.
5. Timer.
6. Absorbent paper

**Handling/Storage:**

1. Reconstitute or dilute only the specific reagents mentioned in the reagent preparation section, when ready to run the assay.
2. Store all kit components at 4°C to 8°C when not in use and do not expose them to temperatures greater than 37°C or less than 2°C
3. Do not use kit components after the expiration date.
4. Do not repeatedly freeze/thaw the reagents as loss of activity may result.

5. Before using, bring all components to Room Temperature (18-25 °C). Upon assay completion return all components to appropriate storage conditions.
6. ELISA plate pouches contain dessicant. Keep the plates sealed in the pouch in the refrigerator when not in use.

**Health Hazard Warnings:**

1. Reagents that contain preservatives may be harmful if ingested, inhaled or absorbed through the skin.
2. Handle Stop Solution carefully. Obtain medical attention in case of accidental ingestion of kit components.
3. Avoid assay of samples containing Sodium azide as it is hazardous.

**Reagent Preparation (all reagents should be diluted immediately prior to use):**

1. 1X Wash Buffer:  
Dilution: To make 1X Wash Buffer, add 50ml of 20X Wash Buffer to 950ml of DI water. This is the working solution.

**Specimen Collection and Preparation:**

Specimens should be clear and non-hemolyzed. Samples should be run at a number of dilutions to ensure accurate quantitation.

**Serum:** Use a serum separator tube and allow clotting for 30 minutes, then centrifuge for 10 minutes at 1000 x *g*. Remove serum layer and assay immediately or store serum samples at < -20°C. Avoid repeated freeze/thaw cycles.

Dilute the serum 1:51 in dilution buffer. For example: add 5 ul of serum sample to 250 ul of dilution buffer. If not assayed immediately, diluted samples should be stored at -20°C or below.

**Procedural Notes:**

1. Read all the instructions thoroughly before performing the test.
2. Allow all reagents to reach Room Temperature before beginning and reconstitute or dilute the required reagents.
3. In order to achieve good assay reproducibility and sensitivity, proper washing of the plates to remove excess unreacted reagents is essential.
4. All Controls and Samples should be assayed at least in duplicates.
5. The assay has been optimized to be used with the protocol mentioned. Any deviation from the same may invalidate the results.

**Assay Procedure:**

1. Bring all reagents to Room Temperature prior to use. It is strongly recommended that all **Controls and Samples** should be run in duplicates or triplicates.
2. Fit the strip holder with the required number of pre-coated Positive Viral Antigen and Negative Viral Antigen. Mark the appropriate strips with a (+) or (-). Allow one well for the Negative Control and one well for Positive Control Sera.
3. Add 50 ul of **Negative Control, Positive Control** and **diluted Sample** to the appropriate wells.
4. Mix the contents in the wells by moving the plate in rapid circular motion, see to it that the contents do not spill. Incubate at **Room Temperature** for **30 minutes**.

5. Aspirate and wash plate 5 times with **Wash Buffer** and blot residual buffer by firmly tapping plate upside down on absorbent paper. Wipe off any liquid from the bottom outside of the microtiter wells as any residue can interfere in the reading step. All the washes should be performed similarly or alternatively a microtiter plate or strip washer may be used.
6. Add **50 ul** of **Anti-Mouse IgG:HRP Conjugate** to the wells.
7. Incubate at **Room Temperature for 30 minutes**.
8. Wash the plate as per the instruction given in step 5.
9. Add **50 ul** of **TMB Substrate** solution into each well of the plate. Incubate for 10 minutes.
10. Add **50 ul** of **Stop Solution** to each well.
11. Measure the optical density of the wells on a plate reader at 450 nm within 15 minutes.

### Interpretation of results

1. It is recommended that each laboratory establish their own criteria for performance of these Research Reagents.
2. In our quality control testing, we use the following criteria:
3. The Positive Control Serum, after subtracting the absorbance in the Negative Control Antigen well, should produce a net absorbance on the Positive Viral Antigen of  $\geq 1.00$  at 450 nm.
4. A sample may be considered Positive by the following criteria: Determine the difference ( $\Delta$ ) 450 nm on the Negative Control Antigen well. This difference ( $\Delta$ ) should be greater than or equal to 1.000 for a sample to be considered Positive.
5. A Value of **1.0 or greater** (after subtracting the absorbance of the Negative Control antigen well) is considered **Positive**. A Value of **less than 1.0** (after subtracting the absorbance of the Negative Control antigen well) is considered **Negative**.

### Expected Values:

The normal value is Negative. Studies have shown that antibodies may take up to 21 days to appear after exposure; therefore, Negative specimen results should be reviewed in relation to a possible exposure date. All Positive specimen results should be confirmed by an alternate method.

### Precautions:

1. Do not mix reagents from different kits or lots. Reagents and/or antibodies from different manufacturers should not be used with this set.
2. Substrate is light and heat sensitive hence do not expose it to direct sunlight while pipetting or incubating.
3. Samples and kit reagents after use should be disposed off observing appropriate regulations.
4. If necessary it is recommended that the results should be confirmed by an alternative method.
5. Do not dilute or adulterate test reagents or use samples not called for in the test procedure.

### LIMITED WARRANTY

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any) from the maker thereof of such non Krishgen made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, and strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

*Krishgen Biosystems. 2023*

***THANK YOU FOR USING KRISHGEN PRODUCT!***

KRISHGEN BIOSYSTEMS®, GENLISA®, DHARMAPLEX™, GENBULK™, GENLISA™, KRISHZYME®, KRISHGEN®, KRIBIOLISA®, KRISHPLEX®, TITANIUM®, QUALICHEK® are registered trademarks of KRISHGEN BIOSYSTEMS. ©KRISHGEN BIOSYSTEMS. ALL RIGHTS RESERVED.

**KRISHGEN BIOSYSTEMS | OUR REAGENTS | YOUR RESEARCH |**